




PRACTICAL INNOVATION

Ziath

DataPaq™ Deep Focus Single Rack Scanner Version 2

Fast, very compact, simple to install and with a 2-year guarantee

The Ziath Deep Focus Single Rack Scanner Version 2 is ideal for tubes that are seated slightly further away from the scanner bed than usual. It can successfully read tubes set on the scanner bed, or up to several millimetres above it.

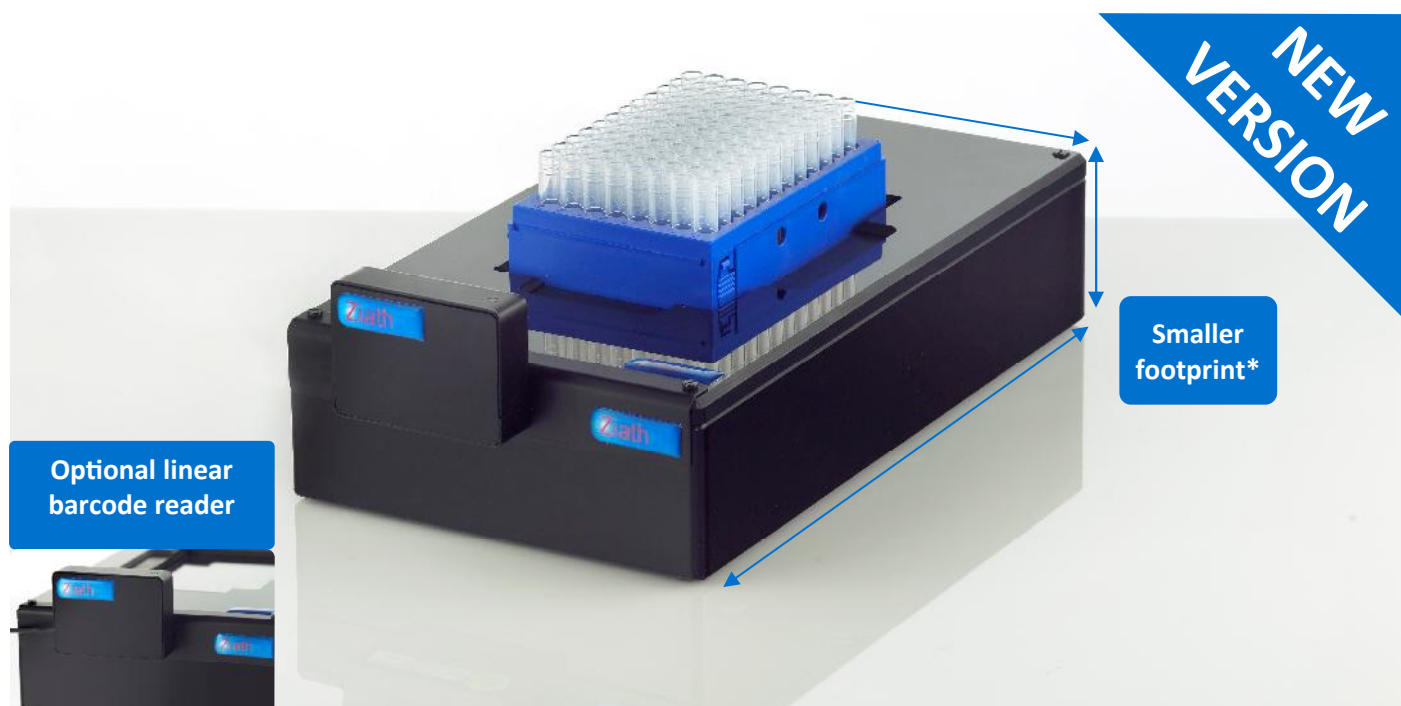

* Smaller than the Ziath DataPaq™ Deep Focus Scanner

Features

- Presence activated - placing rack on scanner activates scan (with linear scanner)
- Compatible with 384, 96, 48 and 24-tube racks
- One-click scan: 10 second scan and decode time per rack
- Integrates fully into robotic platforms
- Compatible with Matrix, ABgene, Nunc, REMP (including upper, locked position), Micronic, FluidX tubes and many more
- Export data to Excel, text, XML, JSON, image
- Operating systems: Windows XP, Vista, 7
- Simple installer; no dongles required
- Database connectivity: SQL Server, Oracle, MySQL, Postgres, HSQL & others
- Very compact: 165 x 288 x 65mm (WxDxH)
- Durable construction - 2 year guarantee
- Free software upgrades for life
- Ziath's expertise and support come as standard



DataPaq™ Deep Focus Single Rack Scanner Version 2

PRACTICAL INNOVATION

Easy Set-Up and Easy To Use

The DataPaq™ Deep Focus Rack Scanner (Version 2) offers easy set-up, straight out of the box, with only a few minutes for configuration.

After a one-time automatic calibration, the scanner is ready to read Matrix, ABgene, Nunc, FluidX, REMP (including upper, locked position) tubes and many other tubes.

Our DataPaq™ software makes it simple to export to Excel, XML, text or JSON and scanned images can also be saved.

The optional Linear Barcode Reader is simply attached to the Deep Focus by a magnet. Once installed, placing a rack on the scanner activates the scan: no mouse-click required.

Intelligent Scan Information

The software displays the results of the scan and can identify the difference between a correctly scanned tube (green) a damaged tube (red) and missing tube (grey square).

Cryoprotection™ Option


The occurrence of condensation whilst scanning tubes straight from cryogenic storage is well known. Ziath's cryoprotective coating eliminates the occurrence and ensures uninterrupted scanning. As Ziath's method for eliminating condensation uses no extra electrical components reliability is increased and costs are significantly decreased.

Ordering Information

Code	Description	Quantity
ZTS-AV62	DataPaq™ Deep Focus Single Rack Scanner Version 2	1
ZTS-AV62-80	DataPaq™ Deep Focus Single Rack Scanner Version 2 with Cryoprotection™	1
ZTS-1DR2-AV6	DataPaq™ Deep Focus Linear Rack Scanner Version 2	1



Ziath's Linear Barcode Reader is easy to install and use.

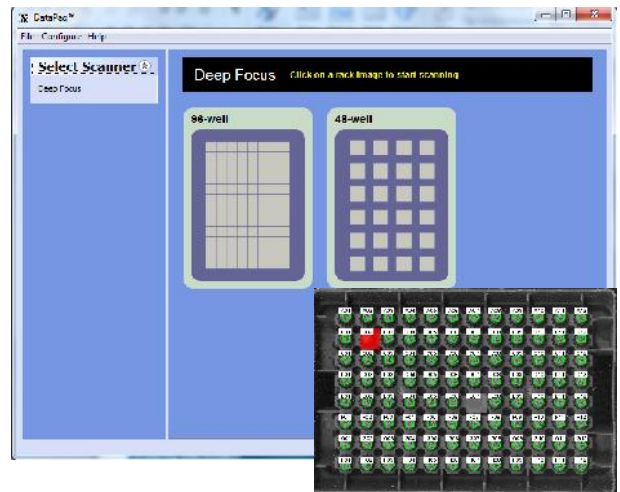

DataPaq™ software: gives intelligent scan information.



Scanner with Cryoprotection™



Scanner without Cryoprotection™